

**BD LSRFortessa H0081
FACSFlow Supply System (FFSS)**

This system automatically supplies sheath from a reservoir in the cart to a pressurized container located on the instrument bench top.

Do not tamper with the pressurized container.

- 1) To refill the *Sheath Reservoir* located on the FFSS cart please do NOT remove the sheath probe or the tank. Fill the reservoir from one of the spare containers that are marked “Sheath” located next to the FFSS. Use the large tank opening for this purpose.**
- 2) To empty the *Waste Reservoir*, carefully remove the probe and place it in the probe holder located on the side of the FFSS next to the waste. Empty the *Waste Reservoir* into the sink and add 100 ml of concentrated bleach. Re-install the container and probe.**
- 3) Leave the FFSS power “ON” at all times. If off due to a power interruption, turn the main switch off and on again.**

Note: If for any reason the FFSS sheath reservoir has been run dry due to customer error, customers must contact UFCR staff to repair the system fluidics.